


# EasyMatch® QC Quick Start Guide

A60-1017-602 ver.1.0

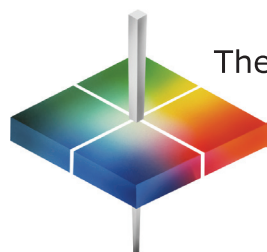

The world's true measure of color

# HunterLab

## A Install EasyMatch QC Software

1. Log into the system using an Administrator account with privileges for the PC — network or local.
2. Insert the installation CD into the CD-ROM drive.
3. Go to Windows **Start > Run > EZMQC\_Menu** and **Open**. The following screen will be shown.

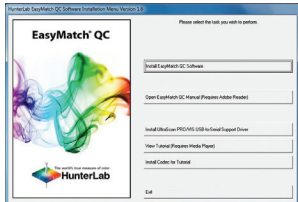


Figure 1. Installation Wizard

4. Select **Install** EasyMatch QC and follow the screen prompts. When prompted, select **SoftKey License** as the type of key to use with the software.

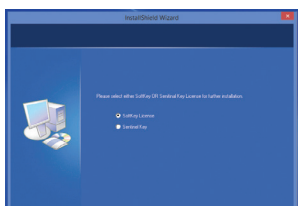


Figure 2. Select SoftKey License

5. When the EasyMatch QC installation is finished, select the button next to **Yes, I want to restart my computer now** and then **Finish** to restart the computer and log back in.

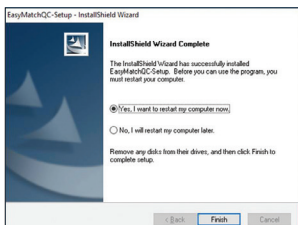


Figure 3. Installation Complete

6. The CD can now be removed.

## B Activate the SoftKey License

1. Select the **EasyMatch QC Icon** or from the Windows Start menu, choose the following to open EasyMatch QC:  
**Start > Programs > HunterLab > EasyMatch QC**

2. A warning message to activate the license will be displayed as shown in Figure 4. Note: EZMQC functions are unavailable before key activation.

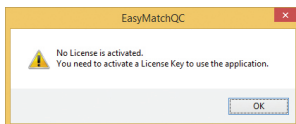


Figure 4. Warning for License Activation

3. The SoftKey License is uniquely associated with the sensor serial number and is provided on a thumb drive supplied with EasyMatch QC or via email from HunterLab.
  - Go to **Help > License Registration/Activation**.
  - Select **Activate License**.

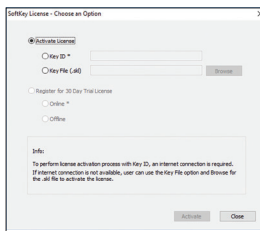


Figure 5. Activate License

### Option #1: Key ID

- a. Copy the License Activation ID from the email or write down the 32-digit code.
- b. From the Choose an Option page (Figure 5), select **Key ID**.
- c. Paste-in or type-in the License Key ID and click **Activate**.
- d. An acknowledgement will be displayed showing the activation status.

Note: An internet connection is required to perform license activation process with Key ID.

### Option #2: Key File

- This method is for using the **Softkey License** (.skl file) on the thumb drive.
- a. Place the thumb drive with the SoftKey License in the USB port.
  - b. From the License Registration page (Figure 5), select **Key File** (.skl).
  - c. **Browse** the USB to find the **SoftKey License** (.skl) file, then click **Activate**.
  - d. An acknowledgement will be displayed showing the activation status.

## C Restart EasyMatch QC

1. When the acknowledgement page has been shown, press **OK** to restart EasyMatch QC.

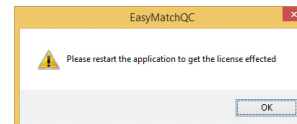


Figure 6. Restart EasyMatch QC

## D Install the HunterLab Sensor.

1. Upon re-start, the following message will be displayed: **'Sensor not yet installed'**. Connect the sensor to the computer using the cable provided. The Setup Sensor screen allows selection of the instrument model and the communications port. To continue select **Add Sensor**. Select **OK** when ready.

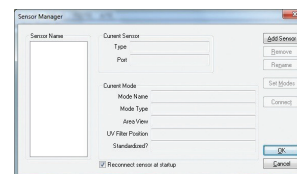


Figure 7. Sensor Manager

2. The next screen sets up the communication port. Select **Next** when ready.

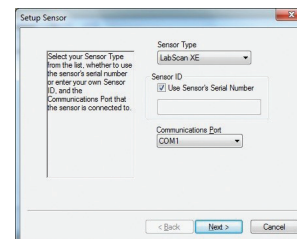


Figure 8. Type & Port Selection

3. Note: The sensor must match the SoftKey License. If a sensor mismatch error occurs, please connect to the correct sensor or contact HunterLab to obtain a license to match the sensor to be connected.

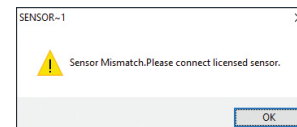


Figure 9. Sensor Mismatch

## F Standardize

1. Select the option next to **Standardize Now** to proceed immediately to standardization upon completion of sensor configuration.

For more detailed instructions on using the instrument, please refer to the EasyMatch® QC User's Manual.

Visit [www.hunterlab.com](http://www.hunterlab.com) for information on all our products.