FAQ: "The slot in the sample clamp on our LabScan XE has worn through. How can we fix this?



We would recommend an upgrade kit that comes with a new spring and sleeve:

A37-1010-689 LSXE Sample Clamp Spring

Used on: LSXE

Spring used in LabScan XE sample clamp.

D04-1010-440 LSXE Sample Clamp Sleeve for Spring

Used on: LabScan XE

Black anodized sleeve used to house LSXE sample clamp spring

You can easily retro-fit this upgrade in the field but unscrewing the little Phillips screw on the underside of the clamp support and re-assembling with the new parts.